

## CELL PELLET PREPARATION FOR RPPA

### Attached Cells:

1. Collect cells by either Trypsin EDTA or scraping cells with rubber police- (minimum of 1 million cells per sample is required)
2. Wash cell pellet twice with PBS (re-suspend and centrifuge)
3. Aspirate PBS as much as possible without disturbing cell pellet
4. Store dry cell pellet in -80 until sample submission
5. *\*\*If shipping samples, pack on plenty of dry ice*

### Suspension Cells:

1. Collect cells by spinning down by centrifugation- (minimum of 1 million cells per sample is required)
2. Wash cell pellet twice with PBS (re-suspend and centrifuge)
3. Aspirate PBS as much as possible without disturbing cell pellet
4. Store dry cell pellet in -80 until sample submission
5. *\*\*If shipping samples, pack on plenty of dry ice*